ecotek 北京力高泰科技有限公司

LI-6800 系统更新操作步骤

(以1.4.02版本为例)

1、首先在电脑上安装 LI-6800-updater-1.4.02 应用程序,下图为软件图标;



- 2、将LI-6800 主机与分析器连接并开机,使用网线连接仪器和电脑;
- 3、双击软件图标 2, 打开软件,单击仪器序列号,再单击 Update LI-6800,弹出升级界面;

LI-6800 Updater (ver 1.4.02) File	- 🗆 X				
Name Model Version 68C-831593 LI-6860 1.3.19 Other	1	🖸 Update LI-6800: 68C-8	31593		1
		LI-6800 Mead (68M-691593)	Version 1.4.0	Needs Update	
		LI-6800 Fluorometer (MPF-631699)	Version 1.4.0	Needs Update	
HAE: 66C-631593 HOST: 66C-631593 IP: 169.254.211.10:6660 IF: fe00::s6d:d80:b8b:716f%18:6660 company: LI-COR Biosciences macaddress: 00:10:94:04:4a:b1 model: LI-6660 sm: 66C-631593 type: Photosynthesis System Console ver: 1.3.16 xver: 1.3.16	2 Update LI-6800	LI-6800 Console (680-631593)	Version 1.4.02	Needs Update	
	What's New in 1.4 Advanced	LI-6800 Soil Chamber (- NA -)	Version 1.6 3	Not Detected	
			Update	Æ	
	Close				

4、单击界面中的 Update, 仪器开始进行升级, 等待 3 个进度条走完, 完成升级。

0 Update U-6800: 680-831	502	2 X	0 Update U-6800: 68C-83	21502	2	
	555	1 ~		1555	1	~
LI-6800 Head (68 H-8 91593)	Version 1.4.0	TRANSFER PROGRESS 2.7%	LI-6800 Head (68H-891593)	Version 1.4.0	SUCCESS!	
LI-6800 Fluorometer (MPF-831699)	Version 1.4.0	TRANSFER PROGRESS 4.4%	LI-6800 Fluorometer (MPF-831699)	Version 1.4.0	SUCCESS!	
LI-6800 Console (68C-831593)	Version 1.4.02	TRANSFER PROGRESS 3.7%	LI-6800 Console (68C-831593)	Version 1.4.02	UPDATING: 01:03	
LI-6800 Soil Chamber (- NA -)	Version 1.6	Not Detected	LI-6800 Soil Chamber (- NA -)	Version 1.6	Not Detected	
	Update			Update		